

PCR-WT Workstation

TOOLS REQUIRED FOR ASSEMBLY

Electric Hand Drill, Phillips Head Bit, 3/8" Magnetic Hex Driver Bit, Socket Set and Ratchet.

BEFORE ASSEMBLING THE WORKBENCH

Read through the assembly instructions to familiarize yourself with the order in which the parts are assembled.

We suggest that you clear & vacuum the area where the bench is to be assembled.

It is important that the bench is assembled in the same sequence as instructed.

HARDWARE

All assembly hardware is provided.




STEP 1

Attach the legs to the modesty panel using (4) 1/4-20 x 1/2" Flange head bolts. The bolts will thread into pre-installed threaded inserts in the sides of the legs




1/4-20 x 1/2"
Flange Head Bolt

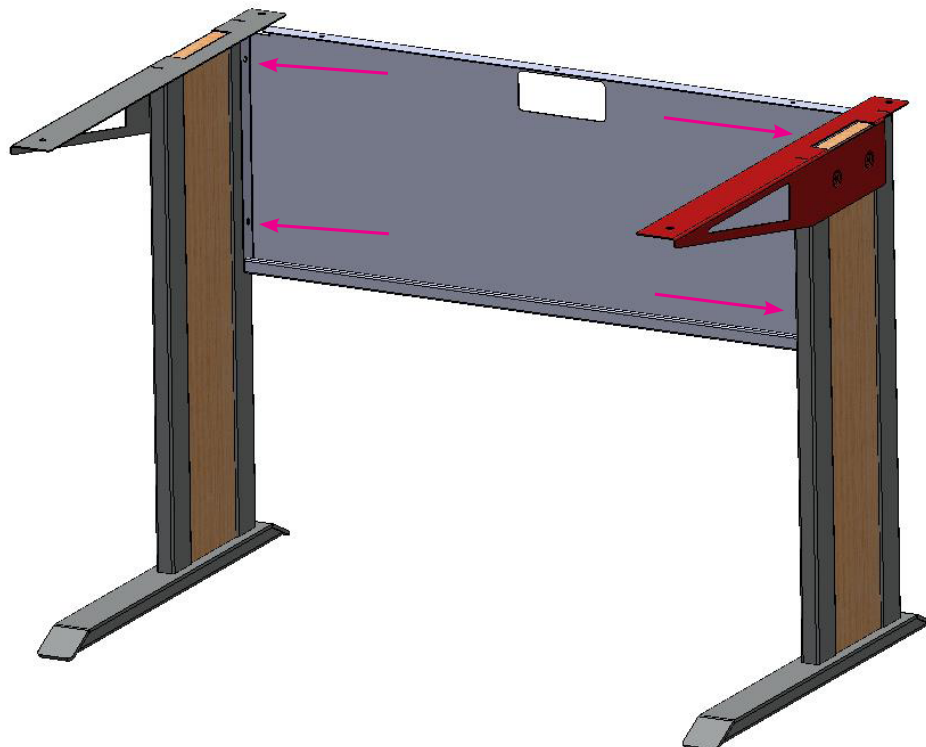

STEP 2

Once legs and modesty panel are attached hang the wire tray over the top of the modesty panel. Holes in the top of the modesty panel will line up with holes in the wire tray.

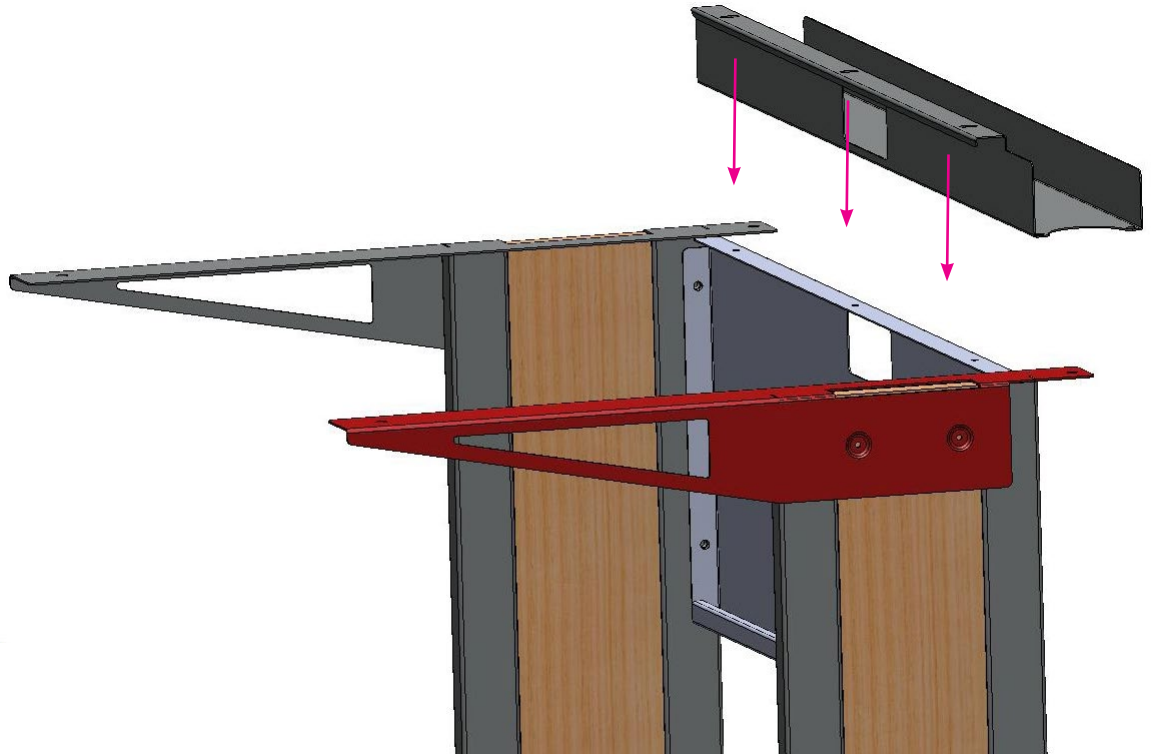

STEP 3

Attach the top to the base. Line up the holes in the legs with the pre-installed wood nuts on the underside of the top. Indicated by pink arrows. Attach using 1/4-20 Flange head bolts. Using Spax wood screws attach the modesty panel/wire tray by screwing through the holes on the top edge. Indicated by green arrows.



1/4-20 x 3/4"
Flange Head Bol



#10 x 5/8"
SPAX Wood Screw

